

SOP Name: T4 Ligase

Author: Marcia Pryce

Source(s): Adapted from Exeter 2015

1. Please note that protocols with Q5 High-Fidelity DNA Polymerase may differ from protocols with other polymerases. Conditions recommended below should be used for optimal performance.

Reaction Setup:

We recommend assembling all reaction components on ice and quickly transferring the reactions to a thermocycler preheated to the denaturation temperature (98°C). All components should be mixed prior to use. Q5 High-Fidelity DNA Polymerase may be diluted in 1X Q5 Reaction Buffer just prior to use in order to reduce pipetting errors.

Materials:

COMPONENT	25 µl REACTION	50 µl REACTION	FINAL CONCENTRATION
5X Q5 Reaction Buffer	5 µl	10 µl	1X
10 mM dNTPs	0.5 µl	1 µl	200 µM
10 µM Forward Primer	1.25 µl	2.5 µl	0.5 µM
10 µM Reverse Primer	1.25 µl	2.5 µl	0.5 µM
Template DNA	variable	variable	< 1,000 ng
Q5 High-Fidelity DNA Polymerase	0.25 µl	0.5 µl	0.02 U/µl

COMPONENT	25 μ l REACTION	50 μ l REACTION	FINAL CONCENTRATION
5X Q5 High GC Enhancer (optional)	(5 μ l)	(10 μ l)	(1X)
Nuclease-Free Water	to 25 μ l	to 50 μ l	

2. Notes: Gently mix the reaction. Collect all liquid to the bottom of the tube by a quick spin if necessary. Overlay the sample with mineral oil if using a PCR machine without a heated lid.

Transfer PCR tubes to a PCR machine and begin thermocycling.

Thermocycling Conditions for a Routine PCR:

STEP	TEMP	TIME
Initial Denaturation	98°C	30 seconds
25–35 Cycles	98°C *50–72°C 72°C	5–10 seconds 10–30 seconds 20–30 seconds/kb
Final Extension	72°C	2 minutes
Hold	4–10°C	